

















DONNING PERSONAL PROTECTIVE EQUIPMENT

While wearing PPE, avoid touching your face or adjusting the PPE once in the infectious area

1	Wash hands	
2	Change into scrubs (if required)	
3	Put on booties (if required)	
4	Put on a lab coat or gown	
5	Put on a head cover (if required)	
6	Put on respirator/ face mask (if required) <ul style="list-style-type: none"> • N-95 respirator for aerosol hazard (must be fit tested annually) • Surgical mask to prevent contamination of experiment 	
7	Put on eye protection or face shield (if required)	
8	Put on gloves – use extended cuff length if available. If the second pair of glove is required, extended pair is preferred for the base layer	

DOFFING PERSONAL PROTECTIVE EQUIPMENT

Remove PPE from the most contaminated to the least

1	Remove booties (if wearing)	
2	Remove gloves <ul style="list-style-type: none"> If wearing double gloves, remove outer glove only 	
3	Remove the lab coat or gown	
4	Wash hands or remove the second pair of gloves	
5	Remove eye protection or face shield (if wearing)	
6	Remove respirator/face mask <ul style="list-style-type: none"> Handle strap only, do not touch the front portion Make sure the mask do not snap back to the face 	
7	Remove head cover (if wearing)	
8	Wash hands	
9.	Change to street clothes (if wearing scrubs) and wash hands	